

#### U-CyTech BV

Yalelaan 48
3584 CM Utrecht
The Netherlands
P +31.30.253 5960
F +31.30.253 9344
INFO@ucytech.com
www.ucytech.com

# Data sheet Rat IFN-y ELISA antibody pair; 20-plate format

Cat. No.: CT700-20

## Coating antibodies (4 vials)

Product: Monoclonal antibody to rat interferon gamma (IFN-y)

Isotype: Mouse IgG<sub>1</sub>

Production: In vitro using serum free medium

Purification: Ammonium sulphate precipitation and DEAE ion-exchange chromatography

Buffer: Prior to lyophilization: 0.25 ml PBS + 125 mM trehalose

Application: Inject 0.25 ml distilled water into the vial and dilute 100 times in PBS. The content

of one vial is sufficient for five 96-well ELISA plates (480 determinations; 50

μl/well).

## Detection antibodies (4 vials)

Product: Biotinylated monoclonal antibody to rat interferon gamma (IFN-γ)

Isotype: Mouse IgG<sub>1</sub>

Production: In vitro using serum free medium

Purification: Ammonium sulphate precipitation and DEAE ion-exchange chromatography

Labeling: With Biotin-7-NHS (N-hydroxysuccinimide)

Buffer: Prior to lyophilization: 0.5 ml PBS + 1% BSA + 125 mM trehalose

Application: Inject 0.5 ml distilled water into the vial and dilute 100 times in PBS + 0.5% BSA +

0.05% Tween-20. The content of one vial is sufficient for five 96-well ELISA plates

(480 determinations; 100  $\mu$ l/well).

#### Standards (10 vials)

Product: CHO-derived recombinant rat interferon gamma (IFN-y)

Application: Cytokine standard in an ELISA system

Reconstitution: Inject 0.5 ml distilled water into the vial. Use immediately.

### Conjugate (4 vials)

Product: SPP conjugate (Streptavidin-HRP)

Application: Inject 0.5 ml distilled water into the vial and dilute 100 times in PBS + 0.5% BSA +

0.05% Tween-20. The content of one vial is sufficient for five 96-well ELISA plates

(480 determinations; 100  $\mu l/well).$ 

The product should be used in combination with TMB substrate.

### General

Sensitivity: 2 pg/ml

Specificity: Validated for detecting natural and recombinant rat IFN- $\gamma$ .

Sterility: Membrane filtered (0.2 µm)

Stability: Lyophilized SPP conjugate is stable for at least one year at -20°C in the dark, the

other lyophilized products are stable for at least one year at 4°C.

After reconstitution, the antibodies are stable for one year at  $4^{\circ}$ C (if kept sterile) and SPP for minimal one year at  $-20^{\circ}$ C in the dark. The reconstituted standard

preparation should be used immediately.

References: Rogóz, Z. et al. 2009. Pharmacol. Rep. 61: 1050-1060

For research use only Version 140822



